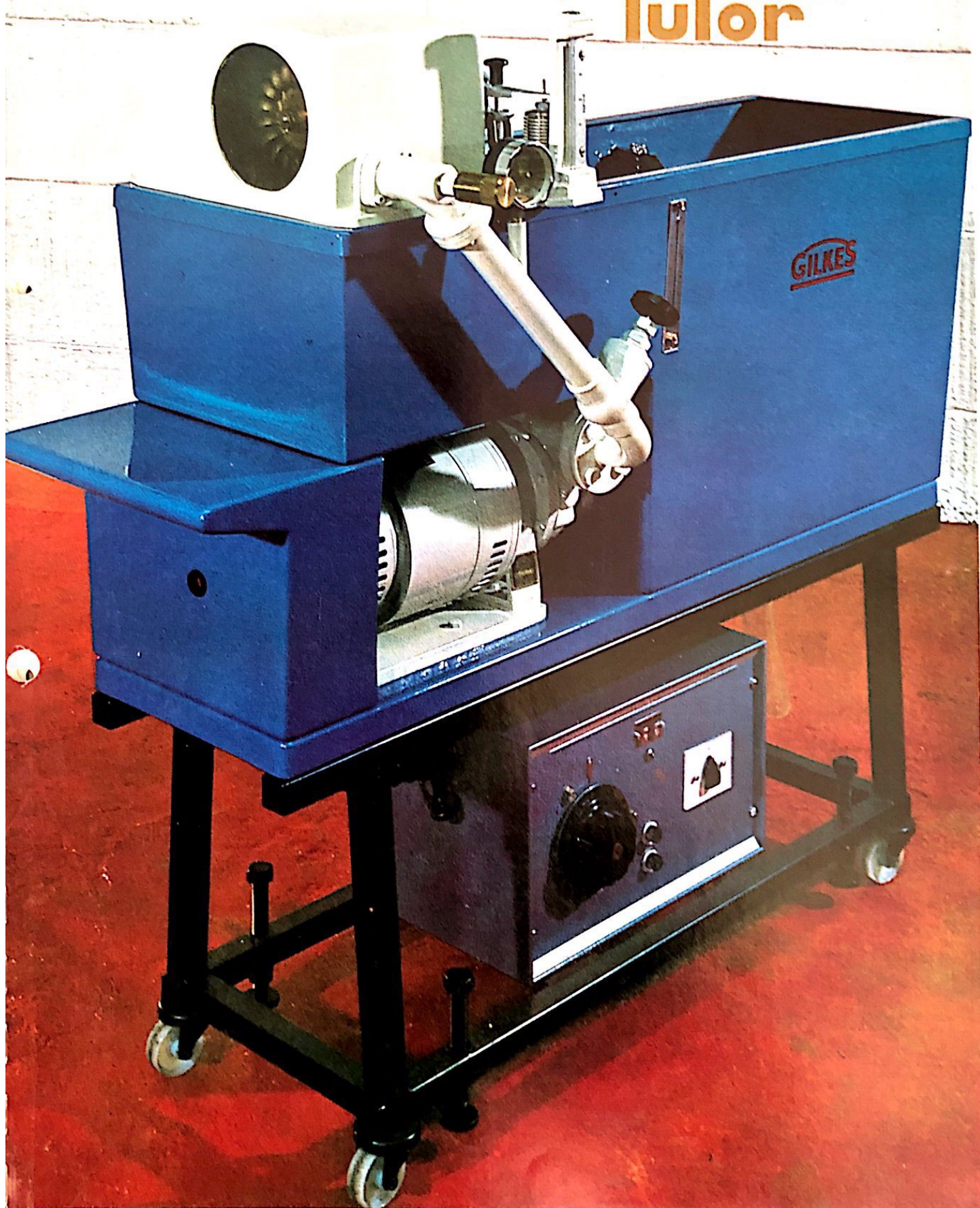
# GILKES Tutor

## INTRODUCTION

The 'Tutor' range of instructional pump and turbine test sets are designed to comply with the trend in modern laboratory design where compactness, adaptability and convenience of handling are important factors. The portability of the units allows storage outside the laboratory or classroom when not in use. 'Tutor' test sets are intended to take their place beside the popular table top fluid mechanics apparatus for flow through orifices, friction loss in pipes, etc. 'Tutor' test sets are of unique value in introducing engineering students to the important concept of 'energy conversion'. Indeed water turbines are the only energy conversion device where the transfer of energy can be seen to take place. It should be noted that a 'Tutor' turbine set fitted with a variable speed pump provides facility for experiments and preparation of iso-efficiency curves for both the pump and the turbine. The results obtained are fully representative of larger pump and turbine units, full details of experiments being given in leaflet M52.

## EACH UNIT COMPRISES:

- **A Pelton wheel or Francis turbine and a motor driven centrifugal pump** which supplies water under pressure to feed the turbine. In the case of the pump set the water is discharged through a throttle valve. **A choice of motors** is available, either **single speed or variable speed**. The variable speed unit incorporates dynamometer mounting. A wide range of supply voltages can be catered for.
- **A fibreglass tank**. The tank comprises two parts: the flow measuring tank and the suction tank. The turbine (or throttle valve) discharges into the upper tank and the water flows over a vee-notch into the lower tank which is directly connected to the pump suction. The rate of flow is measured by means of a water level gauge on the upper tank. The level gauge is calibrated by a scale plate in both English and Metric units. A drain plug is provided.
- **A Prony brake** (for turbine sets only). This is of cast aluminium in two pieces with a heat-resisting brake lining. The brake may be adjusted by means of a handwheel and is particularly smooth in operation. The drum may be cooled with water.
- **A Gilkes force gauge**. This measures the force exerted at the end of the Prony brake arm. A second force gauge is used to measure the load on the dynamometer arm of the motor. In each case the load is applied to the top of a steel bellows. The bellows is filled with mercury and water. As the load is applied mercury is displaced up a glass tube; the height to which the mercury rises indicates the load applied. The force gauges are calibrated in English and Metric units.
- **A pressure gauge**. A dual scale pressure gauge is provided to measure the inlet pressure to the turbine and the discharge pressure from the pump.
- **A trolley with levelling screws**. The trolley makes the 'Tutor' fully portable and supports the test set at a convenient height. The four levelling screws allow maximum accuracy during flow measurements.
- **A motor starter or speed controller**. The single speed version is provided with a direct-on-line starter mounted near the motor. The variable speed version is provided with a transformer/rectifier speed controller. Infinitely variable speed control is possible and the unit will operate from a wide range of AC supply voltages.

## INTERCHANGEABILITY

The Pelton Wheel and Francis turbine units are easily interchangeable. The conversion kits may be purchased separately, if required. If a 'Tutor' pump test set has been purchased it may be converted to a turbine test set by the purchase of a Francis or Pelton conversion kit. Many principals will prefer to purchase two complete sets and so save the time taken to change from one turbine to another.

## SPEED MEASUREMENT

No tachometer is fitted. The speed should be measured with a hand tachometer that will be readily available in any laboratory. Each exposed shaft end has a suitable centre. A hand tachometer can be supplied as an optional extra.

## STROBOSCOPE

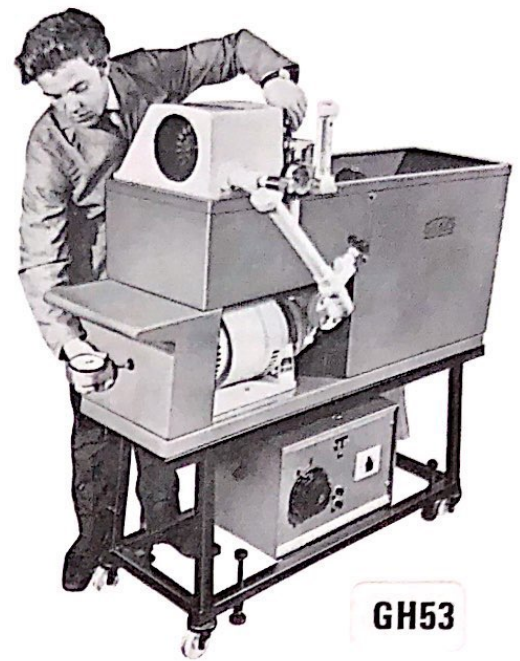
An electronic stroboscope may be supplied as an optional extra. This unit may be used to show the flow patterns in both Francis and Pelton turbines. In addition the unit is calibrated in revolutions per minute and may thus be used to measure the speed of the turbine. Full descriptive literature will be supplied upon request.

## POLYTHENE COVER


A polythene cover is provided with each 'Tutor' set.

## CHARACTERISTIC CURVES


Sets of detailed characteristic curves are available on request.

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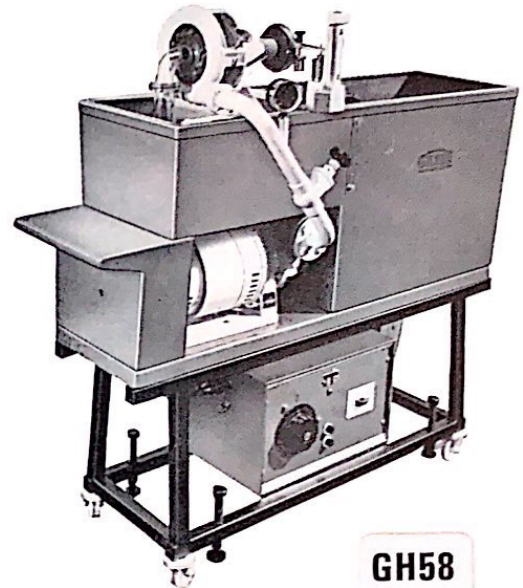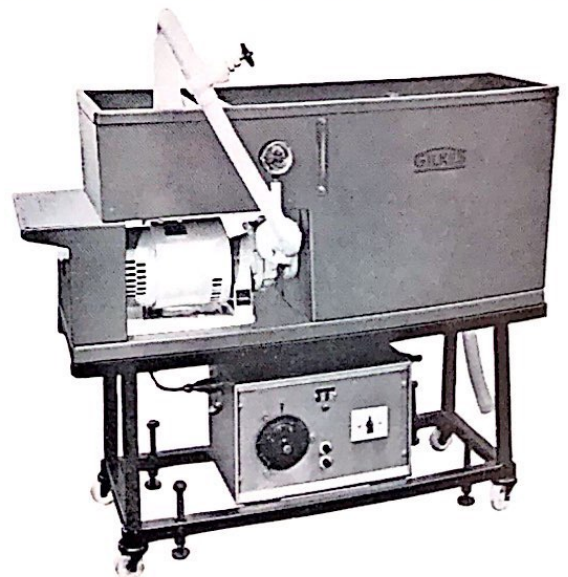# PELTON

(Items 50, 51, 52 & 53)

This small impulse turbine is an exact replica of modern industrial units of the best design. The maximum efficiency is approximately 70%. Experiments may be conducted to draw the students' attention to the similarity which exists between impulse water turbines and impulse steam turbines. Use of a stroboscope vividly demonstrates the water flow patterns on entry into and discharge from the buckets.

## SPECIFICATION

### OUTPUT POWER:

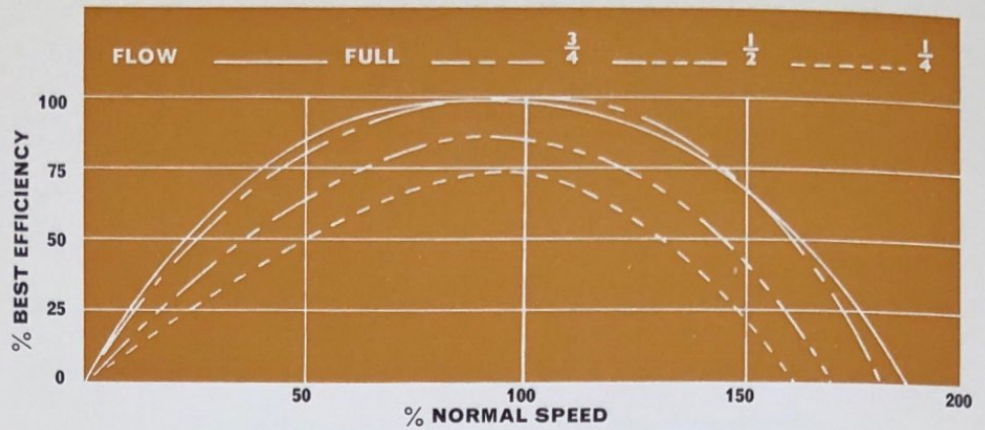
0.3-0.38 k.W. at 1600 r.p.m. on 21 m. head.

### CASING:

Rust proof fibreglass with plexiglass window.

### RUNNER:

Precision bronze casting.

### BEARINGS:

Deep groove ball, sealed to give protection from corrosion.

### SHAFT:

Stainless steel.

### SPEAR & NOZZLE ASSEMBLY:

Gunmetal, with calibrated micrometer hand-wheel to adjust the position of the spear valve. Repeatability of experiments is thus assured.

### SPECIFIC SPEED:

Metric 24.3, English 5.6.

# FRANCIS

(Items 54, 55, 56 & 58)

This small reaction turbine, in spite of its small size, has all the features, hydraulic and mechanical of modern industrial units of the best design. The maximum efficiency is approximately 65%. The guide vanes are adjustable and may be seen through the plexiglass window and draft tube. Experiments may be conducted to draw the students' attention to the similarity between a reaction water turbine and reaction steam turbine.

## SPECIFICATION

### OUTPUT POWER:

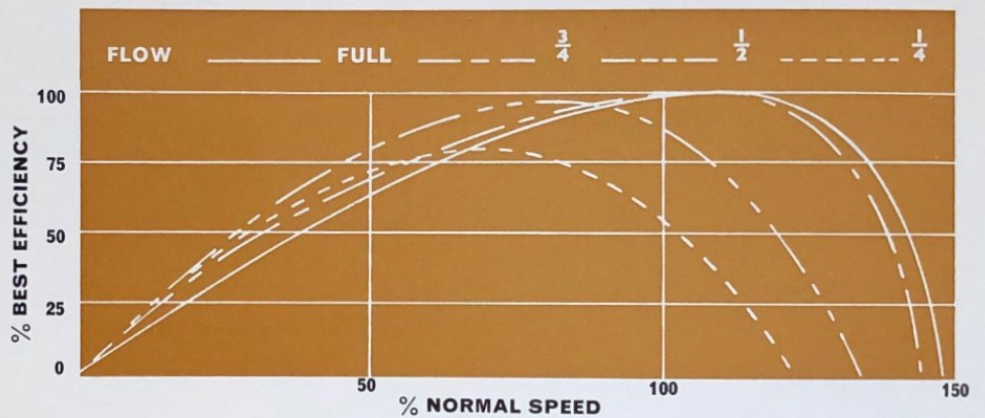
0.3-0.38 k.W. at 2350 r.p.m. on 15 m. head.

### CASING:

Bronze casting.

### RUNNER:

Precision bronze casting.

### BEARINGS:

Deep groove ball, sealed to give protection from corrosion.

### SHAFT:

Stainless steel.

### GUIDE VANES:

Bronze, fully adjustable with calibrated setting scale. Repeatability of experiments is thus assured.

### END COVER & DRAFT TUBE:

Plexiglass assembly.

# PUMP

The centrifugal pump is fitted to all 'Tutor' sets and may be driven by a choice of electric motors. In the case of the 'Tutor' Pump Test Sets (Items 60 & 61) a dynamometer motor is always provided which enables full iso-efficiency curves to be determined. It should be noted that the pump is of generous size, yielding truly representative characteristic curves.

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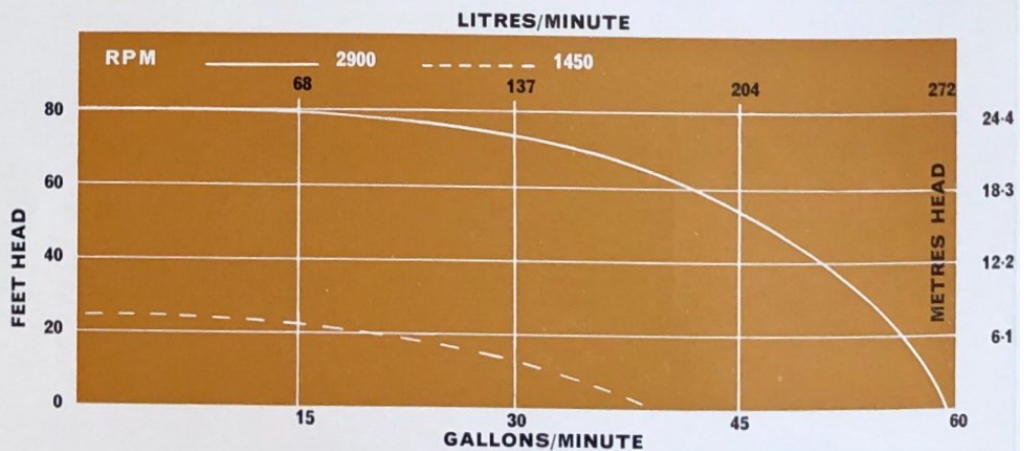
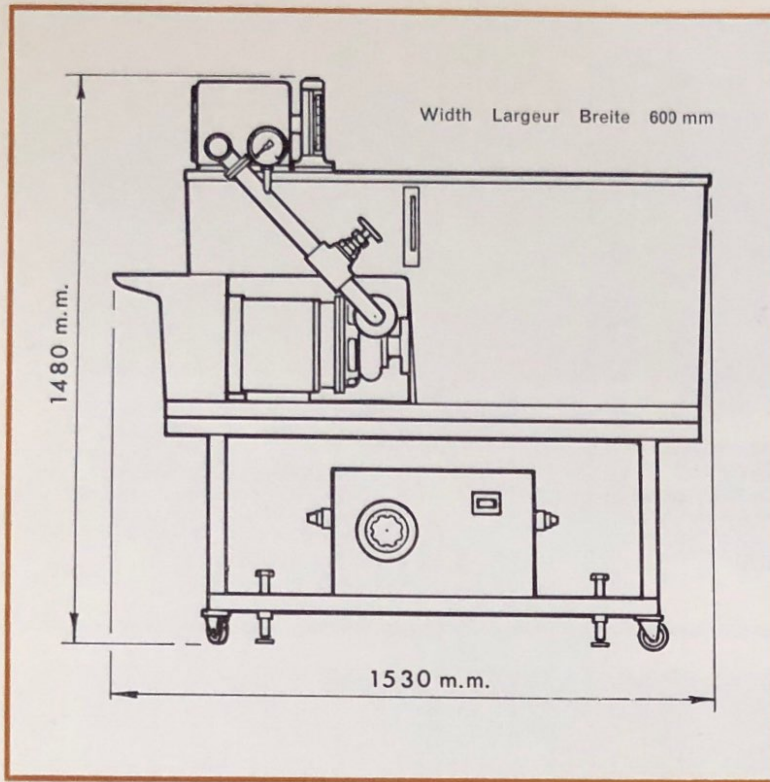
Fig. 4 Dimensions of 'Tutor' Pelton set with variable speed motor (Item 53). All other models are similar in size.

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**APPROXIMATE SHIPPING SPECIFICATION FOR ALL STANDARD UNITS**

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Modifications may be made to design and data to incorporate latest developments. The information given in this leaflet is, therefore, liable to alteration without notice.

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**Other products in the Gilkes range include**

- Larger Pelton turbines
- Francis pump-turbines
- Axial flow pump-turbines
- Flow visualisation channel
- Parallel-series pump set
- Kaplan turbines
- Gas turbine test set
- Special centrifugal pump

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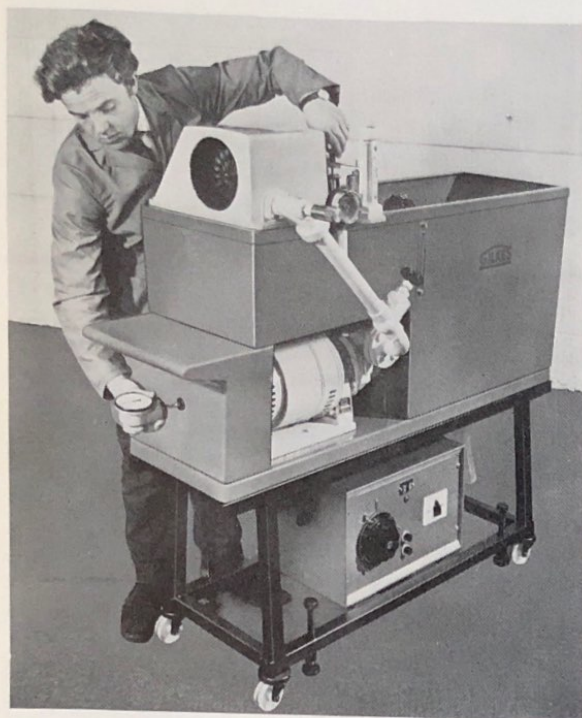
Fig. 1. Tutor Pelton Unit with variable speed motor. (Item No. 53).

### PELTON TURBINES

Two Pelton turbines of different sizes are available – the 'Tutor' and the 'No. 12'. Together with accessories these units offer a wide choice of specifications including fully self-contained closed-circuit systems. Larger Pelton turbines with governing facilities are available for special applications.

#### 'Tutor' Pelton

The Gilkes 'Tutor' Pelton is complementary to the 'Tutor' Francis. The unit is completely self-contained and is supplied complete with supply pump, pipes, 'V' notch weir, instruments, fibreglass tank and trolley. A choice of electric motor drives for the supply pump is available and of these the variable speed version is the most useful since full tests may then be conducted on the centrifugal pump as well as the turbine. An interchangeable Francis turbine unit is available. Construction features include: pump and turbine made entirely of bronze and stainless steel, corrosion-proof fibreglass tank and a plexiglass inspection window for the turbine. Despite its modest dimensions this machine can provide valuable test results and curves. For further details ask for leaflets M23 and M52.

#### 'No. 12' Pelton

This is a larger turbine having a 240 mm mean diameter compared with the 100 mm Tutor runner. The turbine can operate on supply heads up to 60 m when a power of 6 kw is developed. The turbine may be supplied alone (Item 1) or with a closed circuit system and supply pump (Item 62). A large plexiglass window is provided permitting observation of the runner and water jet. Standard instrumentation consists of prony brake, pressure gauge, force gauge and tachometer.

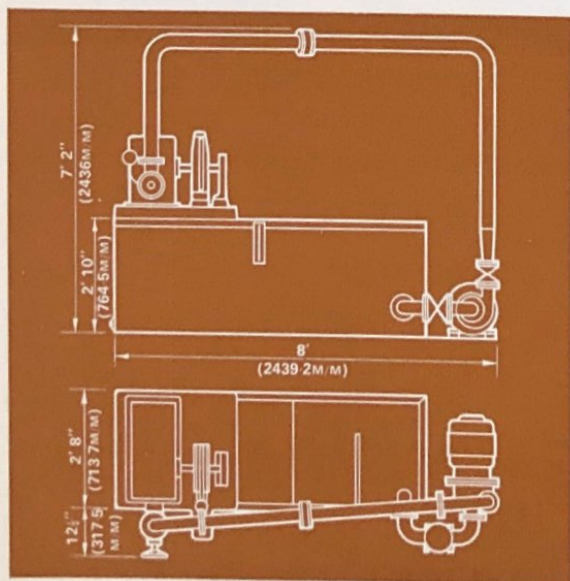
Fig. 2. No. 12 Pelton closed circuit system. (Item No. 62).

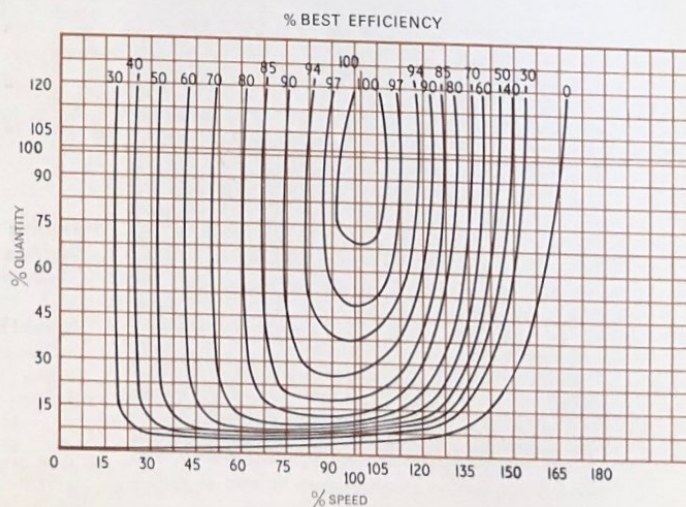
Fig. 3. Pelton Iso Curve Efficiency/Speed Flow.

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## TURGO IMPULSE TURBINES

This design of medium head impulse turbine is made only by Gilkes. The jet of this turbine is set at an angle to the face of the runner and is discharged at the opposite side. With a specific speed of 11.5 rpm English units or 51.5 rpm metric units, it extends the advantages of the impulse turbine into the lower specific speed portion of the Francis turbine range. The Turgo Impulse Wheel is now used in many important installations throughout the world. It can be supplied with a complete closed circuit system.

1 JET/SPEAR TRAVEL. THEORETICAL. 3 % EFFY/%Q No. 12 PELTON  
2 JET/SPEAR TRAVEL. ACTUAL. 4 % EFFY/%Q TUTOR PELTON

Fig. 8. Impulse Turbine and nozzle characteristics.

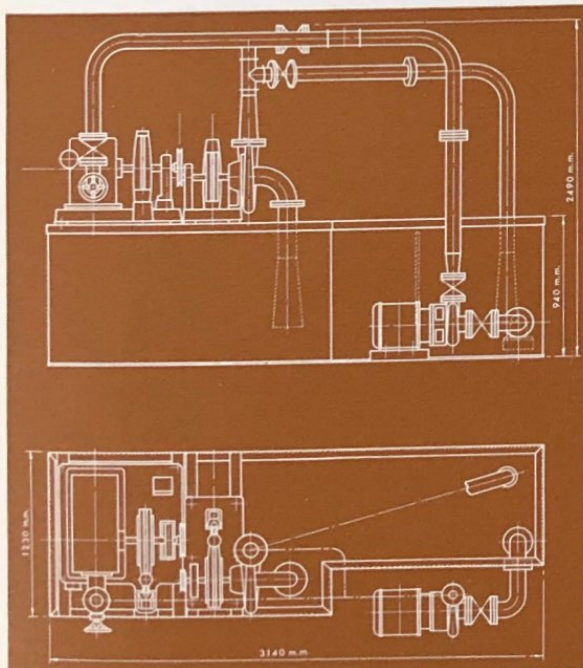
Fig. 7. Closed circuit system incorporating Francis & Pelton Turbines. (Item No. 64).

	Tutor	No. 12		7.5" T.I.	
		Std.	Max.	Std.	Max.
Net head feet	70	50	200	50	100
Net head metres	21.4	15.3	61	15.3	30.5
Flow: igpm	34	84	169	326	462
litre/sec.	2.5	6.4	12.7	24.6	34.9
Normal speed	1600	650	1290	766	1085
Runaway speed	2550	1015	1435	1380	1955
Power: bhp	0.4/0.5	1.03	8.0	3.8	10.8
Specific speed:					
English	5.6	5.1		11.3	
Metric	24.3	22.6		50.2	
(L/S & Metres)					

Fig. 9. Impulse Turbine data.

**GILKES**

**GILBERT GILKES & GORDON LTD.**

KENDAL · ENGLAND · Telephone: KENDAL 20028 · Telex: 65125

Printed in England 2M/11 72 1969

# EDUCATIONAL WATER TURBINES

## Francis Type

GILBERT GILKES & GORDON LTD

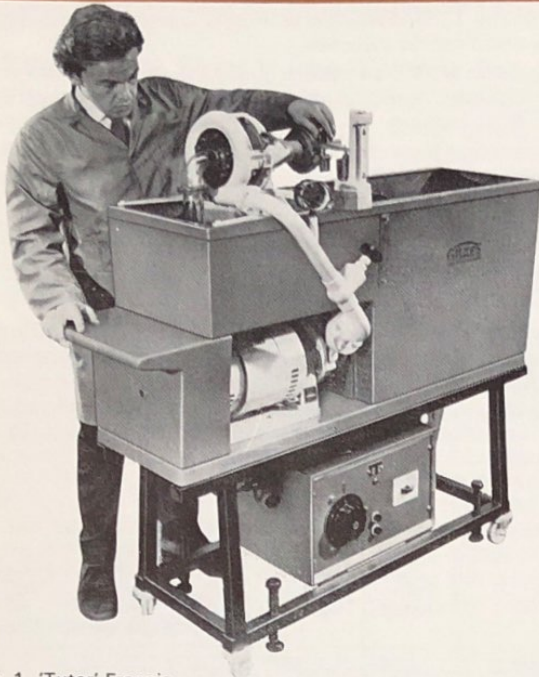
Fig. 1. 'Tutor' Francis unit with variable speed dynamometer motor. (Item No. 58).


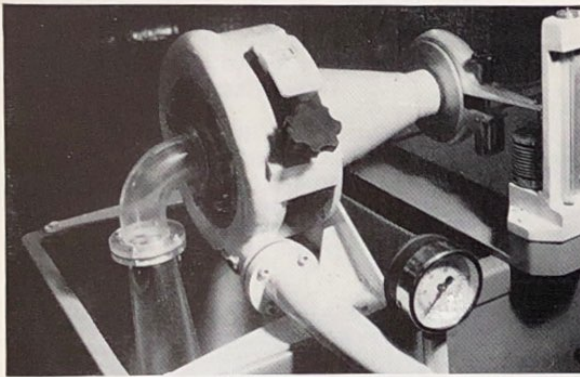
Fig. 2. Detail of 'Tutor' Francis turbine showing guide vane adjusting mechanism.
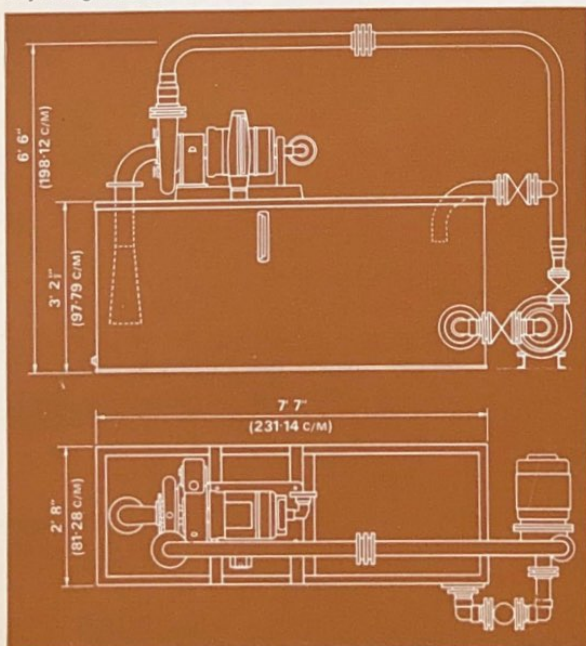
### FRANCIS TURBINES

Two Francis turbines are available: the 'Tutor' and the 150 mm R.P.T. (Reversible Pump Turbine). The latter machine has a runner of 150 mm diameter and is of advanced specification which will, in addition to normal Francis turbine experiments provide curves of performance whilst operated as a centrifugal pump. The machine is available in several versions and is described in greater detail overleaf.

### 'Tutor' Francis

The Gilkes 'Tutor' Francis is complementary to the 'Tutor' Pelton. The unit is completely self-contained and is supplied complete with supply pump, pipes, 'V' notch weir, instruments, fibreglass tank and trolley. A choice of electric motor drives for the supply pump is available and of these the variable speed version is the most useful since full tests may then be conducted on the centrifugal pump as well as the turbine. An interchangeable Pelton turbine unit is available. Construction features include: pump and turbine made entirely of bronze and stainless steel, corrosion-proof fibreglass tank and a plexiglass inspection window for the turbine. Despite its modest dimensions this machine can provide valuable test results and curves. For further details ask for leaflets M23 and M52.

### 150 mm Reversible Francis Pump Turbine

Designed as a modern pump-turbine this unit has many special features. Three principal versions are available:—

- With friction (Prony) brake. (Item 16).
- With dynamometer motor-generator. (Item 18).
- With dynamometer motor-generator and hydraulic governor (Item 19).

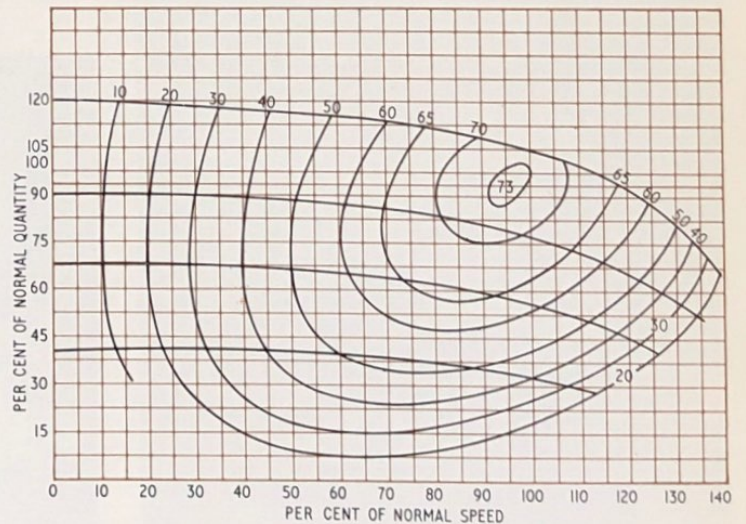
Fig. 4. Iso-efficiency curve for 'Tutor' Francis turbine.

← Fig. 3. 150 mm Reversible Francis pump/turbine on tank, fitted with motor generator and governor. (Item No. 67).

*Our policy is one of continuous improvement and we reserve the right to alter any details of our products at any time without giving notice.*

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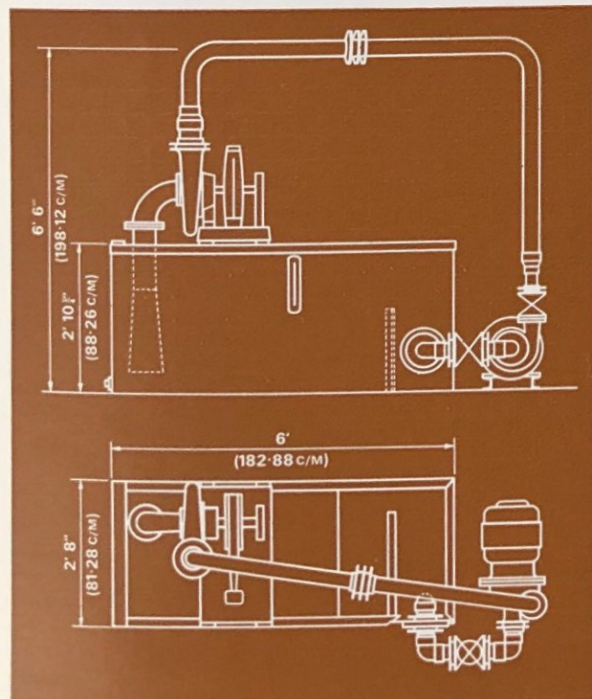
Fig. 7. 150 mm 'R.P.T.' on tank with pump. (Item No. 63).

As an extra, each of the three versions may be supplied with a plexiglass inspection window and open impeller. Pressure tappings are provided in the plexiglass and energy transfer across the impeller may thus be studied. Refer to Items 16P, 18P and 19P. More technical details of these turbines are available in leaflet M48.

The wide range of experiments includes:—

- Proof of the hydrodynamic similarity between centrifugal pumps and Francis turbines.
- Investigation of the 'complete quadrant' of the machine as a normal pump, normal turbine, reverse rotation pump and reverse flow turbine.
- Characteristics during governor control of guide vanes.
- Typical Francis turbine characteristics.
- Typical centrifugal pump characteristics.
- The effect of variable diffuser vanes in a centrifugal pump.

### Closed Circuit Systems

We strongly recommend that turbines be installed in closed circuit form as shown in Figs. 3 and 7 since this permits maximum flexibility in the laboratory and guarantees satisfactory operation. The systems are supplied complete in every respect and require only simple assembly, filling with water and connection to an electricity supply for immediate operation. (Refer to Items 63, 66 and 67).

### DUTIES

	'Tutor' Francis	Turbine 150 mm 'RPT'	150 mm 'RPT' with Motor Genr.	150 mm 'RPT' with Motor Genr. Govr.	150 mm 'RPT' as Pump
Net Head : feet	50	70	70	70	52
metres	15.3	21.4	21.4	21.4	16
Flow : igpm	45	218	218	218	149
litre/min	204	980	980	980	660
Normal speed : rpm	2360	2000	2000	2000	2200
Runaway speed : rpm	3850	3500	3500	3500	—
Power : bhp	0.4/0.5	3.4	3.4	3.4	3.28
Specific speed :					
English	12.6	18.2	18.2	18.2	1387
Metric, $Q = m^3/sec.$	54.7	80.8	80.8	80.8	29.4
Gross weight : lb/kg	532/240	450/210	952/440	1030/470	850/390
Volume : cu.ft/cu.m	29/0.82	28/0.80	28/0.80	34/0.97	28/0.80

Fig. 8. Technical data for instructional Francis turbines.

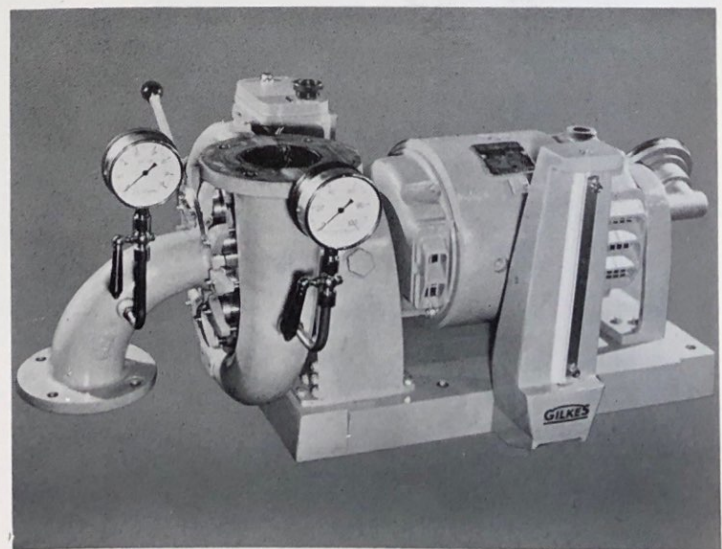
Fig. 9. 150 mm 'R.P.T.' Reversible pump/turbine fitted with motor generator, oil pressure governor, tachometer and force gauge. (Item No. 19).

**GILKES**

**GILBERT GILKES & GORDON LTD.**

KENDAL · ENGLAND · Telephone: KENDAL 20028 · Telex: 65125

Printed in England 2M/11/72